

## Chemical compatibility of CHROMAFIL® products

The chemical compatibility depends on several parameters such as time, pressure, temperature and concentration. In most cases, CHROMAFIL® filters will have only short contact with a solvent. In these cases they may be used despite of limited compatibility. For example, a PTFE filter with PP housing does not liberate any UV-detectable substances during filtration of 5 mL THF, although PP shows only limited resistance towards THF. The following table lists the chemical compatibility of our CHROMAFIL® materials.

Solvent	Material											
	MV	CA	RC	PA	PTFE	H-PTFE	PVDF	PES	PET	GF	IC	PP
Acetaldehyde	-	-	+	O	+	+	+	+	+	+	+	O
Acetic acid, 100 %	-	-	-	-	+	+	+	+	+	+	+	+
Acetone	-	-	+	+	+	+	-	-	+	+	+	+
Acetonitrile	-	-	+	+	+	+	+	+	+	+	+	+
Ammonia, 25 %	-	-	O	-	+	+	+	+	O	+	-	+
Benzene	+	+	+	+	+	+	O	+	+	+	+	O
n-Butanol	+	+	+	O	+	+	+	+	+	+	+	+
Cyclohexane	+	+	+	O	+	+	+	+	+	+	+	+
Dichloromethane	+	-	+	-	+	+	+	-	+	+	-	-
Diethyl ether	O	O	+	+	+	+	+	+	+	+	+	O
Dimethylformamide	-	-	O	+	+	+	-	-	+	+	+	+
1,4-Dioxane	-	-	+	+	+	+	O	-	+	+	+	O
Ethanol	-	+	+	+	+	+	+	+	+	+	+	+
Ethyl acetate	-	-	+	+	+	+	+	+	+	+	+	O
Ethylene glycol	O	O	+	+	+	+	+	+	+	+	+	+
Formic acid, 100 %	+	-	O	-	+	+	+	+	O	+	+	+
Hydrochloric acid, 30 %	-	-	-	-	+	+	+	-	+	-	-	+
Methanol	-	-	+	+	+	+	+	+	+	+	+	+
Nitric acid, 65 %	-	-	-	-	O	+	O	O	O	+	-	-
Oxalic acid, 10 % aqueous	+	-	+	-	+	+	+	+	+	+	+	+
Petroleum ether	+	+	+	+	+	+	+	+	+	+	+	+
Phosphoric acid, 80 %	-	-	O	-	+	+	O	+	+	+	-	+
Potassium hydroxide, 1 mol/L	-	-	O	+	+	+	O	O	O	+	+	+
2-Propanol	+	+	+	+	+	+	+	+	+	+	+	+
Sodium hydroxide, 1 mol/L	-	-	O	+	+	+	O	O	O	O	+	+
Tetrachloromethane	+	-	+	+	+	+	O	+	+	+	+	O
Tetrahydrofuran	-	-	+	O	+	+	+	-	+	+	+	O
Toluene	+	-	+	+	+	+	+	+	+	+	+	O
Trichloroethene	+	+	+	O	+	+	+	O	+	+	+	O
Trichloromethane (chloroform)	+	-	+	-	+	+	+	-	+	+	+	-
Urea	+	+	+	+	+	+	+	+	+	+	+	+
Water	+	+	+	+	+	+	+	+	+	+	+	+
Xylene	+	+	+	+	+	+	O	O	+	+	+	O

Data not guaranteed.

+ resistant, - not resistant, O limited resistance

## Material

### Membranes

MV = cellulose mixed esters, CA = cellulose acetate, RC = regenerated cellulose, PA = polyamide,

PTFE = polytetrafluoroethylene, H-PTFE = hydrophilized polytetrafluoroethylene, PVDF = polyvinylidene difluoride, PES = polyethersulfone,

PET = polyester, GF = glass fiber, IC = special filter for ion chromatography

### Housing material

PP = polypropylene

[www.mn-net.com](http://www.mn-net.com)

**MACHEREY-NAGEL**



MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6-8 · 52355 Düren · Germany

DE / International:

Tel.: +49 24 21 969-0

Fax: +49 24 21 969-199

E-mail: info@mn-net.com

CH:

Tel.: +49 62 388 55 00

Fax: +41 62 388 55 05

E-mail: sales-ch@mn-net.com

FR:

Tel.: +33 388 68 22 68

Fax: +33 388 51 76 88

E-mail: sales-fr@mn-net.com

US:

Tel.: +1 484 821 0984

Fax: +1 484 821 1272

E-mail: sales-us@mn-net.com

